

## 1. Product: Spheroid Dish (Cat. No. 111350)


SPL Spheroid Dish is designed to easy culturing and counting the spheroid. Meshes inside the well facilitate identifying and counting the spheroid. After experiment, the spheroid can be easily collected.



## 2. Experimental condition

- ❖ Cell line: A549(Human epithelial cell), HeLa(Human epithelial cell), ADSCs(Human mesenchymal stem cell)
- ❖ Cell Density:  $5 \times 10^5$  cells / 800  $\mu$ l
- ❖ Culture Condition: 37°C, 5% CO<sub>2</sub>

## 3. Results

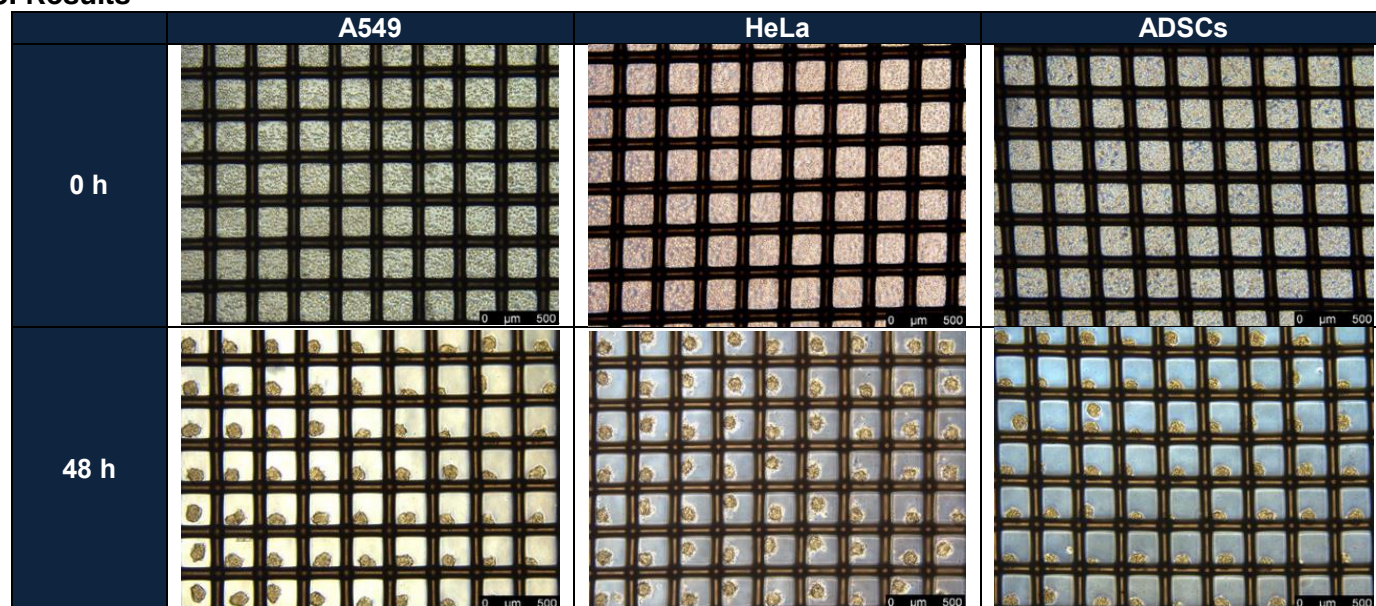

## 4. Conclusion

Spheroid Dish showed the result of cell spheroid formation with high uniformity for various cells.

## 5. Ordering Information

Cat.No.	Material (Dish / Mesh)	Dish Style d x h (mm)	Internal Well Dimension d x h (mm)	Low Binding Treatment	Sterile	Packaging
111350	PS / PE	35.00 x 10.00	25.00 x 1.00	+	+	3 / 15

[www.spllifesciences.com](http://www.spllifesciences.com)

For technical assistance, contact SPL R&D Center at:  
Tel: +82-31-533-4800; Fax: +82-31-533-1430; e-mail: [spl@ispl.co.kr](mailto:spl@ispl.co.kr)

To place an order, contact your local distributor or  
Tel: +82-31-533-4800; Fax: +82-31-533-1430; e-mail: [business@ispl.co.kr](mailto:business@ispl.co.kr)